


Micro-Pak Mold Screening Test: Sampling Procedure



1. Remove the slide from the provided plastic case.

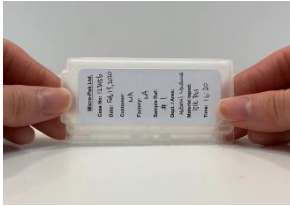

2. Record the sample location, sample details on the slide and label provided, paste the label at the back side of the plastic case.



3. Peel off protective liner from the slide to expose the adhesive.

Note: Do NOT touch the sampling window. The adhesive is very tacky and only needs to be pressed gently. Do NOT press down hard.




4. Place the slide with the adhesive side down on the surface to be sampled. **(Use the centering line to orient the adhesive over the exact sample location desired).** Gently press down to ensure contact is made following the below procedure:

- For samples with **no visible mold, 5 - 6 areas** of one sample should be collected on one tape. Please pay attention to the pressure of each collection, do not over press. (Note: Try to cover all main, lining and different materials of the sample)
- For samples with **visible/suspected mold, up to 3 areas** of one sample should be collected on one tape to avoid over collection.



5. Carefully remove the slide from the surface and place it back into the plastic case **(Sampling window should be facing UP).**

Note: Do NOT replace the plastic "peel" liner.

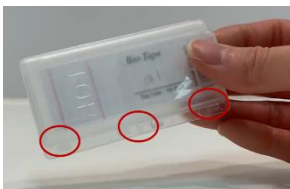

6. Close the slide case properly.

Note: Ensure it is firmly closed on the three circled areas in red.



7. Put the slide case in a Ziploc bag and send it to the nearest Micro-Pak Technical Office together with the provided "Mold Screening Test Application Form" for further analysis.